



Everything you wanted to know about Benchtop NMR (but were afraid to ask)

SNUG 2024 Postgraduate Course in
NMR Spectroscopy

3rd Dec. 2024

Robin J. Blagg PhD MChem MRSC


Applications Team Leader – Magnetic Resonance
Oxford Instruments

Introduction


- + Benchtop NMR
- + Benchtop NMR Spectroscopy
- + Small Molecule Characterisation / Analysis
- + Reaction Monitoring & FlowNMR
- + Time-Domain NMR



What is NMR ?

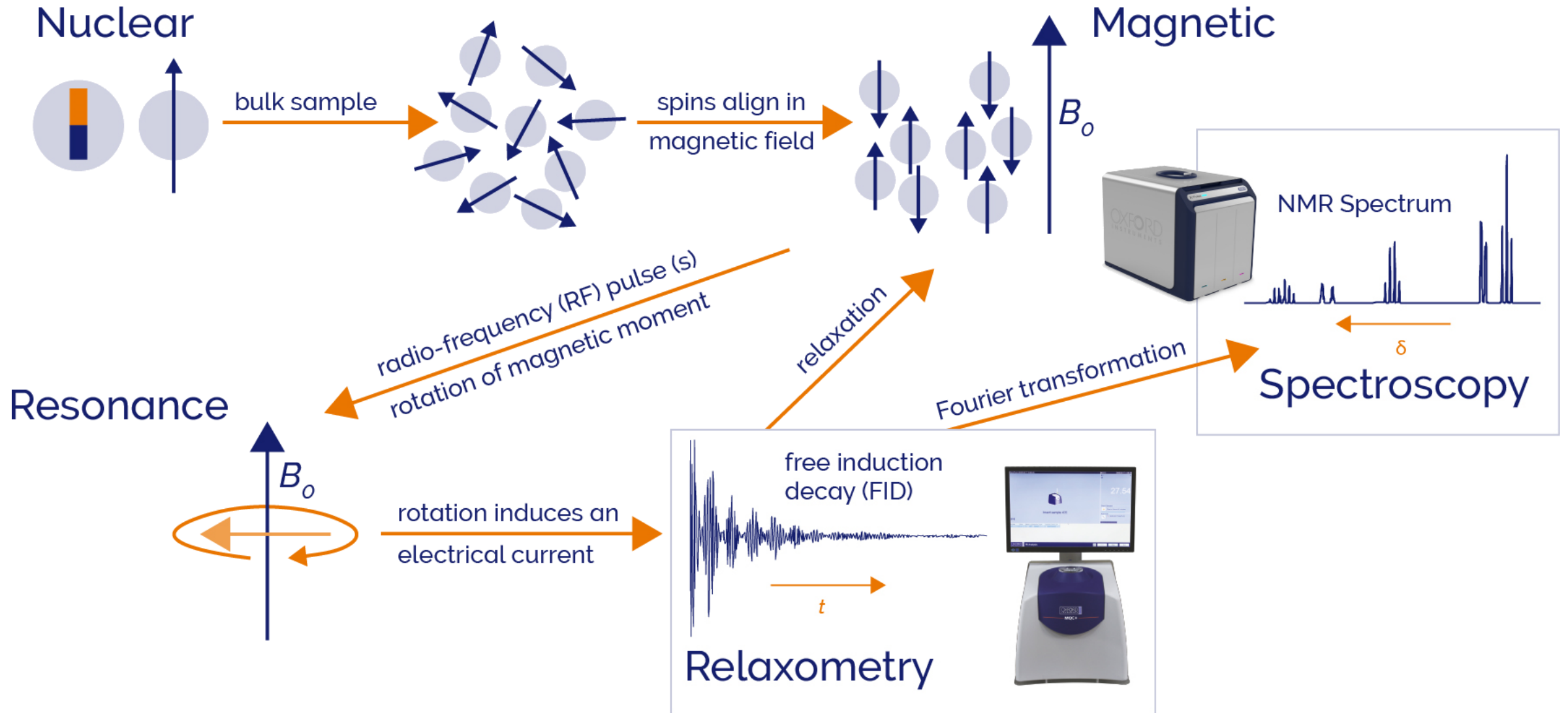


What is an NMR Spectrometer / Relaxometer ?

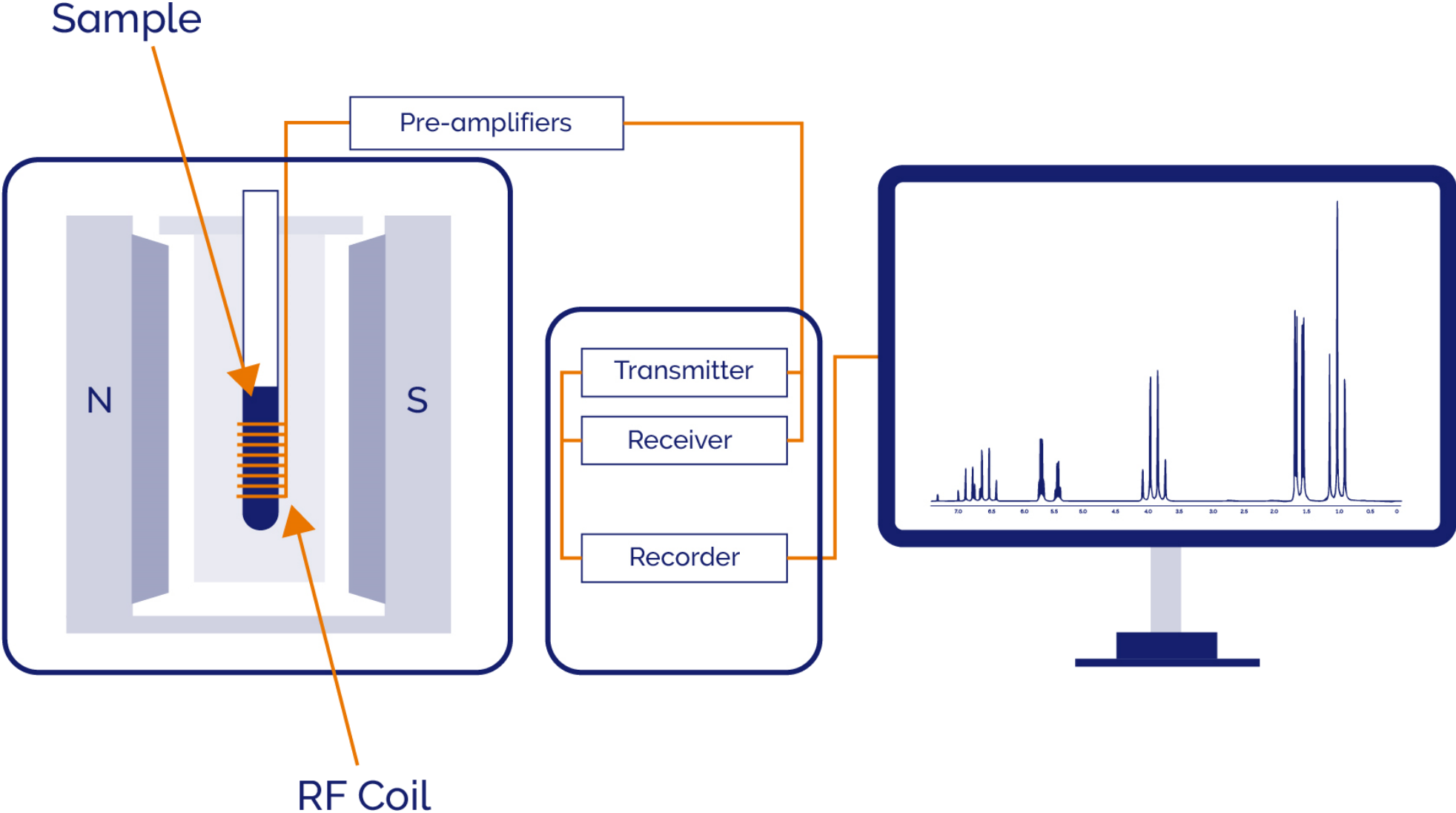


Benchtop versus High-field NMR

- Benchtop NMR
 - up to 120 MHz
 - permanent magnets ($\text{Nd}_2\text{Fe}_{14}\text{B}$)

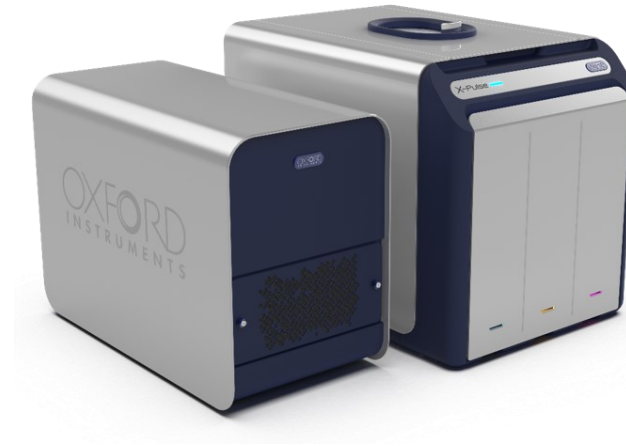

- High-field NMR
 - up to 1.2 GHz (28.2 Tesla)
 - superconducting magnets
 - liquid cryogenes (N_2 & He)
 - dedicated facilities / personnel

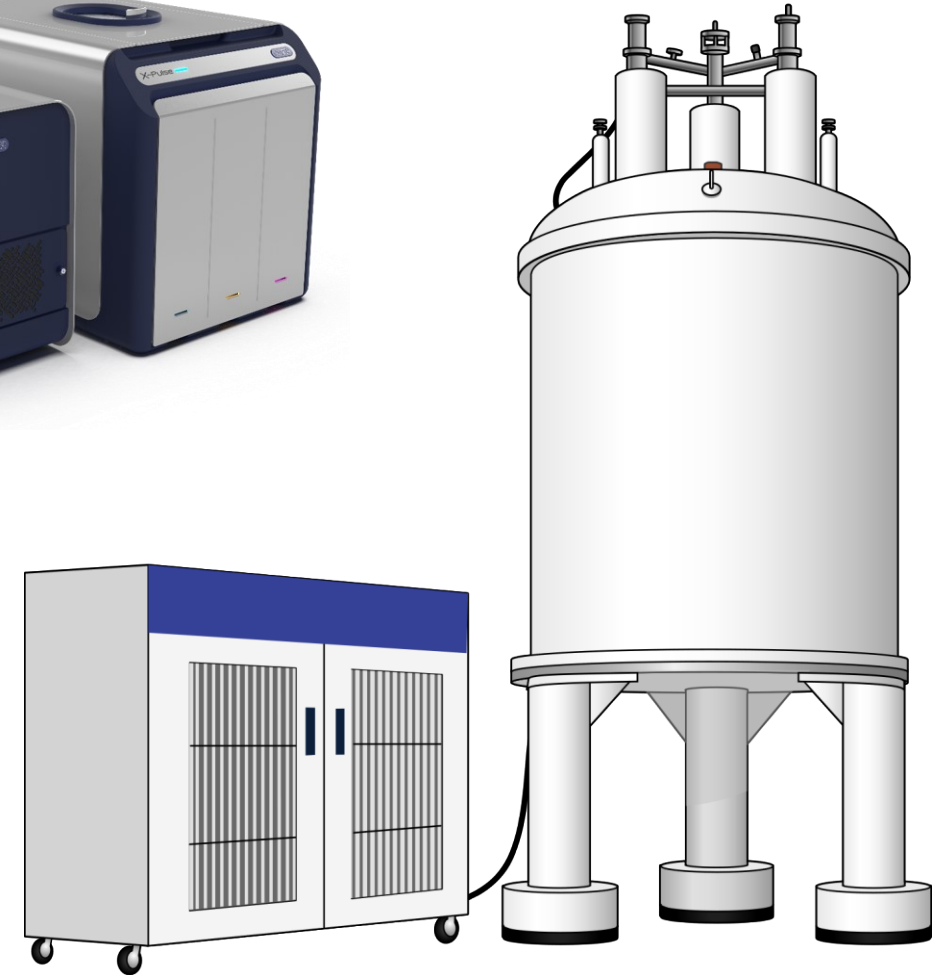

X-Pulse

Broadband Benchtop NMR Spectrometer

X-Pulse Broadband Benchtop NMR Spectrometer




- 1.4 T / 60 MHz permanent magnet
- High resolution (CHCl_3 in acetone- d_6)
 - $\nu_{1/2} \leq 0.35 \text{ Hz} / \leq 10 \text{ Hz} @ ^{13}\text{C}$ satellites
- High sensitivity (1% ethylbenzene in CDCl_3)
 - ≥ 140 signal-to-noise ratio (SNR)
- Three-axis field gradients as standard (z-axis: $\geq 0.5 \text{ T}\cdot\text{m}^{-1}$)
 - gradient-selective two-dimensional homo- & hetero-nuclear correlation
 - pulsed field gradient spin echo, for diffusion measurements
- Shaped pulses and gradients as standard
 - solvent suppression using WET (**W**ater suppression **E**nhanced through **T**₁ effects)
 - selective one-dimensional TOCSY, NOESY

X-Pulse Broadband Benchtop NMR Spectrometer

- Three channel spectrometer
 - Proton/fluorine channel
 - Broadband X-channel (*optional*)
 - ^7Li , ^{11}B , ^{13}C , ^{23}Na , ^{27}Al , ^{29}Si , ^{31}P , and many more ...
 - Deuterium lock channel
- User removable / exchangeable probe
- Auto-sampler, 25 positions (*optional*)
- Variable temperature operation (*optional*)
 - minimum range 0 to +65°C
- Flow NMR capable (*optional*)
- Mobile workstation/cart



- Sequences programmed in Python
 - complete control of spectrometer hardware
 - pre- / post-acquisition processing



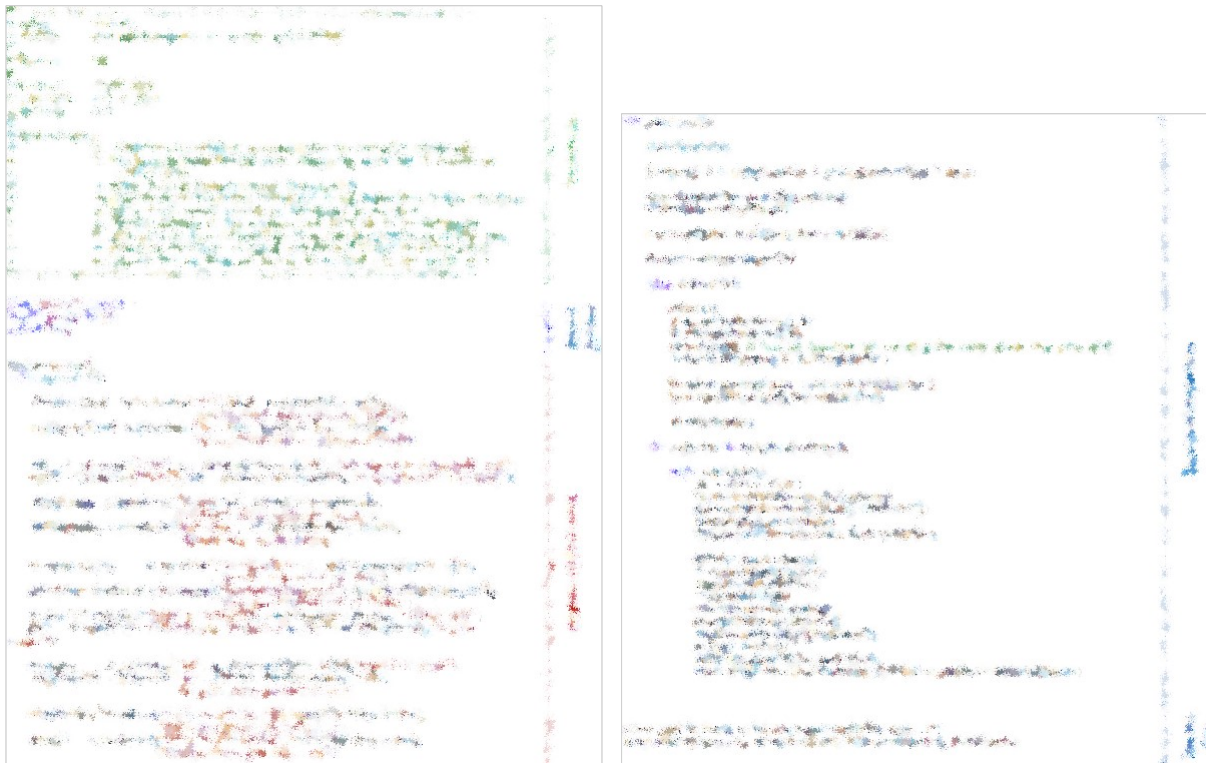
The image shows two side-by-side screenshots of Python code. The left screenshot displays a complex script with various comments and function calls, likely related to spectrometer hardware control. The right screenshot shows a more structured code block with comments in German, such as 'Erzeugung der Pulse', 'Ausgabe der Parameter', and 'Ausgabe der Ergebnisse', indicating the generation and output of pulse sequences.

- Two radio-frequency channels
 - high- & low-power
 - shaped / adiabatic pulses
- Three bipolar gradient channels
- External (TTL) triggers
 - input & output
- Customisable JCAMP-DX output files





Small Molecule Characterisation / Analysis

Small Molecule Characterisation / Analysis





- Characterisation & analysis of pure compounds & mixtures
 - ^1H , ^7Li , ^{11}B , ^{13}C , ^{19}F , ^{23}Na , ^{27}Al , ^{29}Si , ^{31}P , ^{59}Co , ^{195}Pt ...
 - one- & two-dimensional pulse sequences
 - FID, InvRec, CPMG, INEPT, DEPT, PGSE ...
 - COSY, TOCSY, NOESY, HSQC, HMBC, HMQC ...

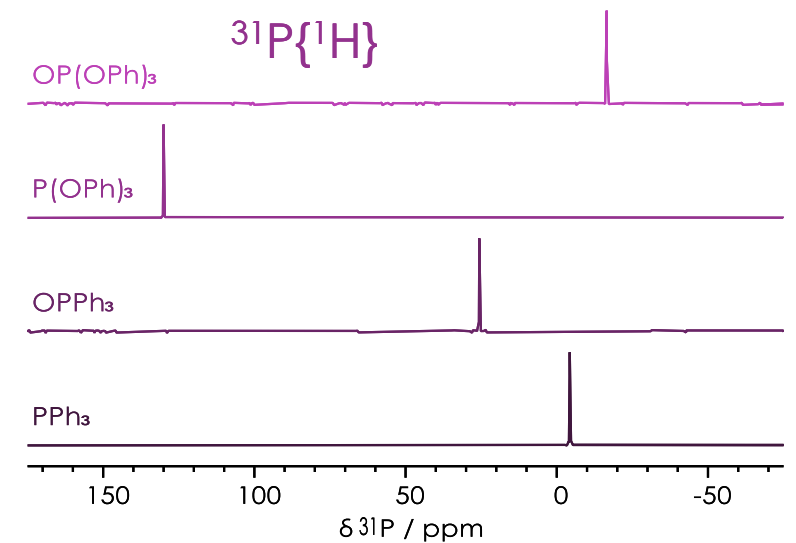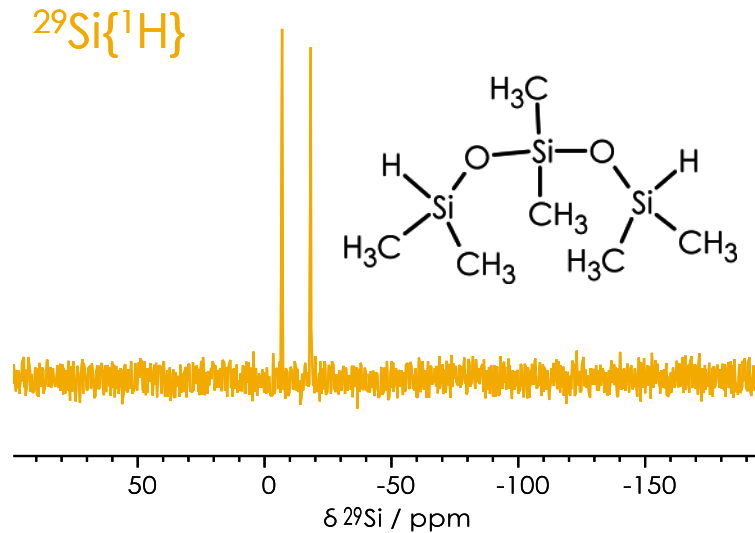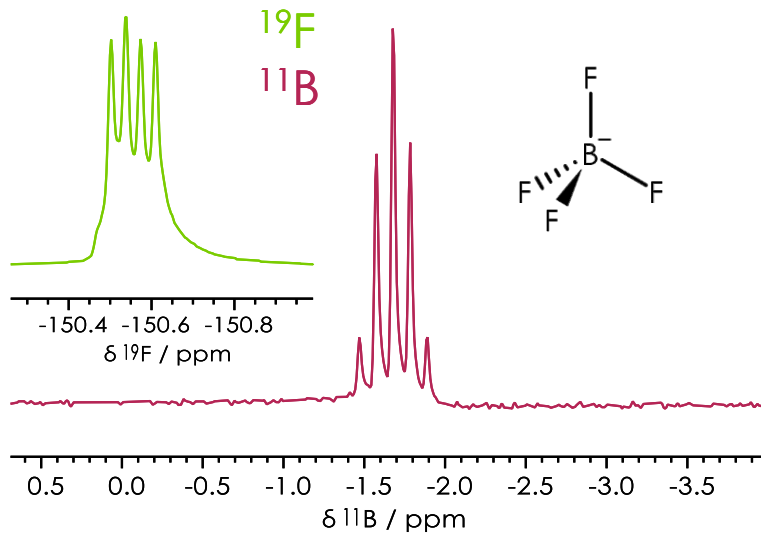





Small Molecule Characterisation / Analysis

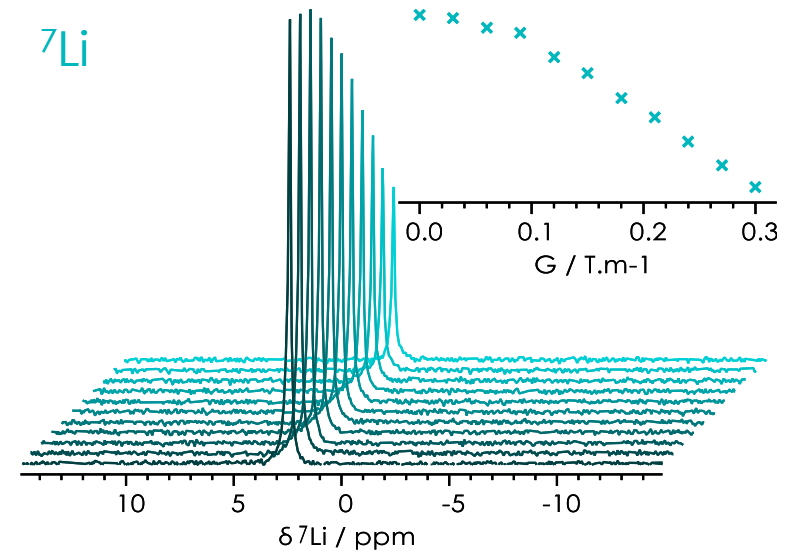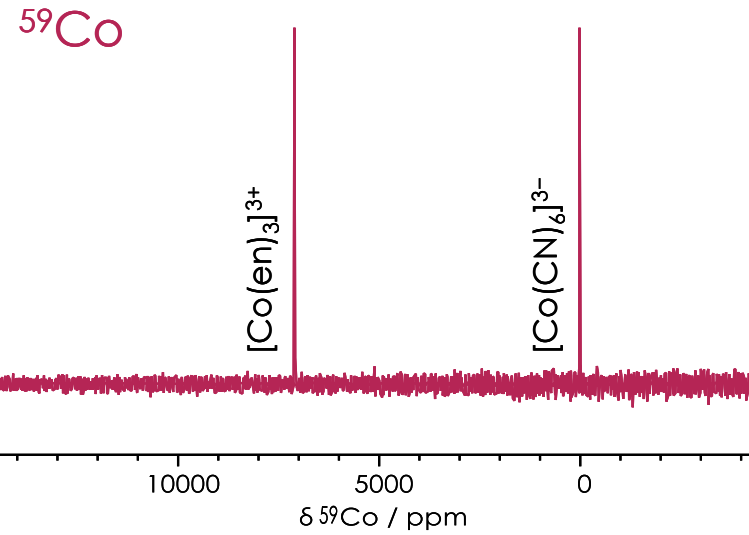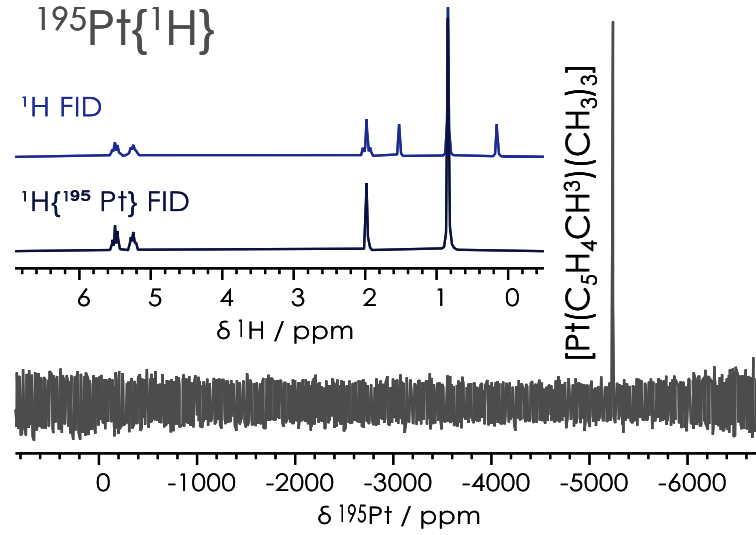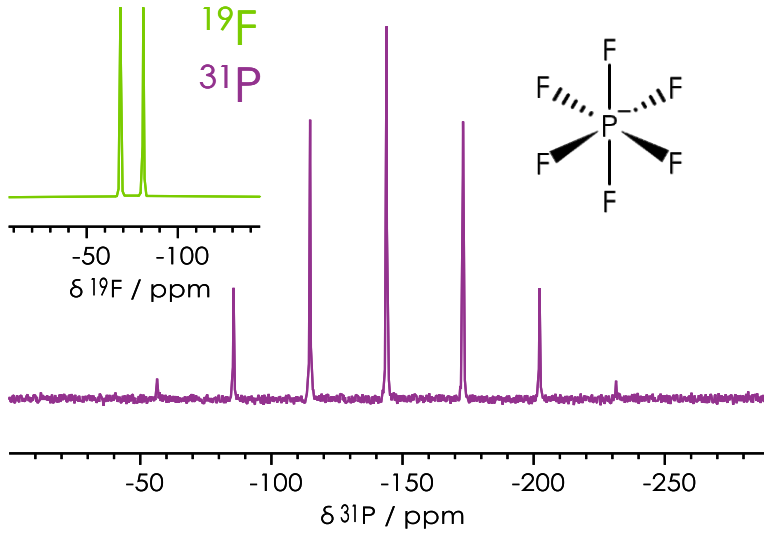

Small Molecule Characterisation / Analysis

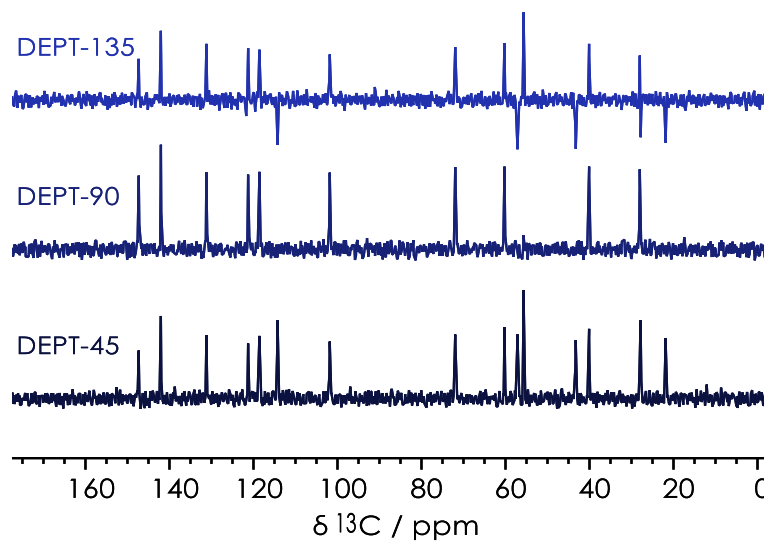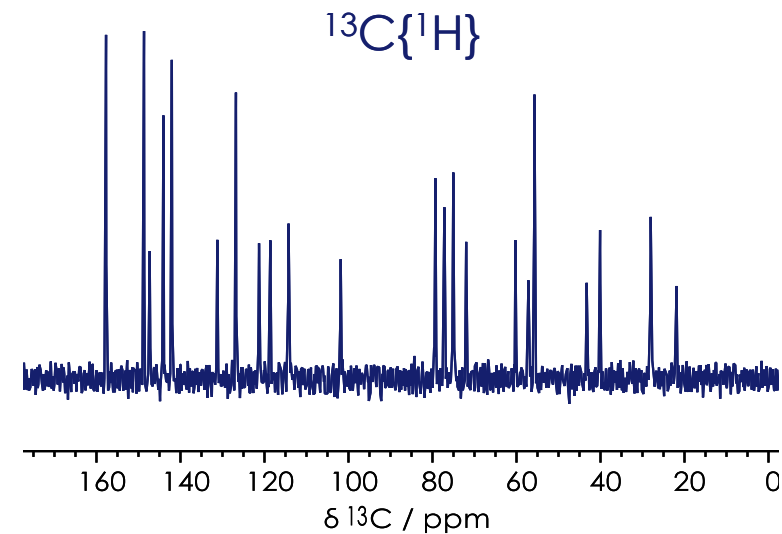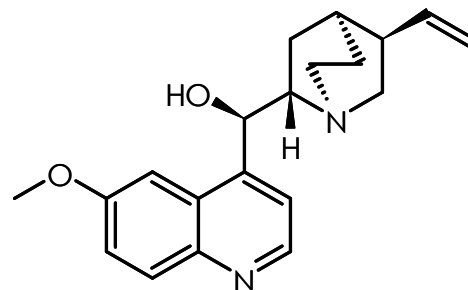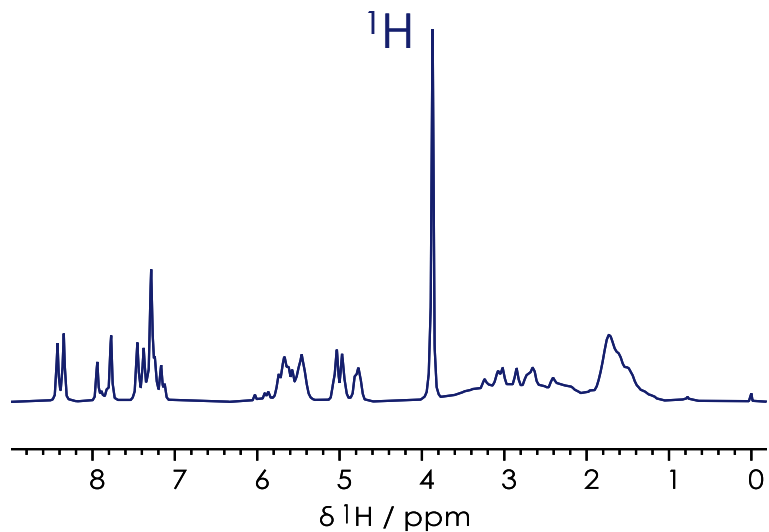


Small Molecule Characterisation / Analysis

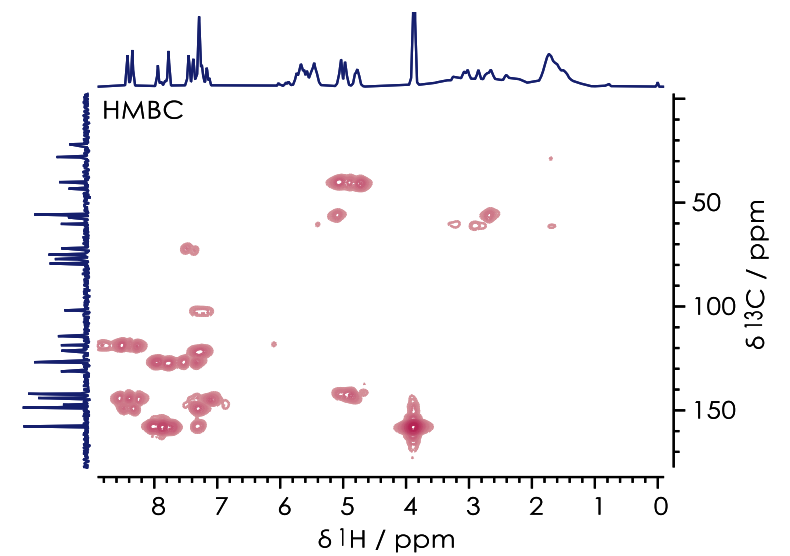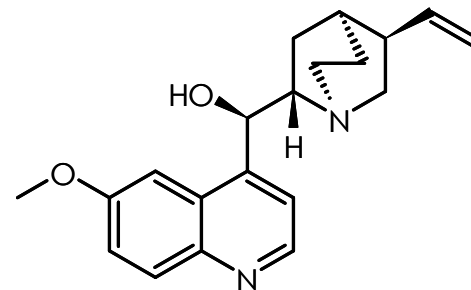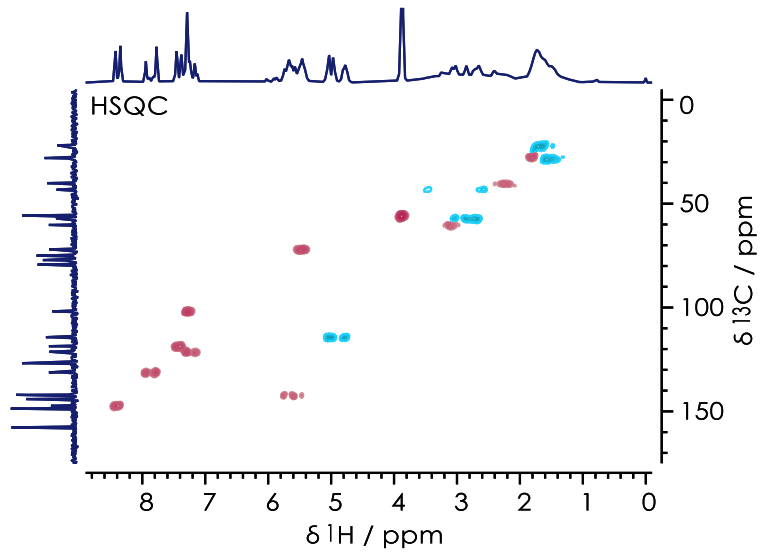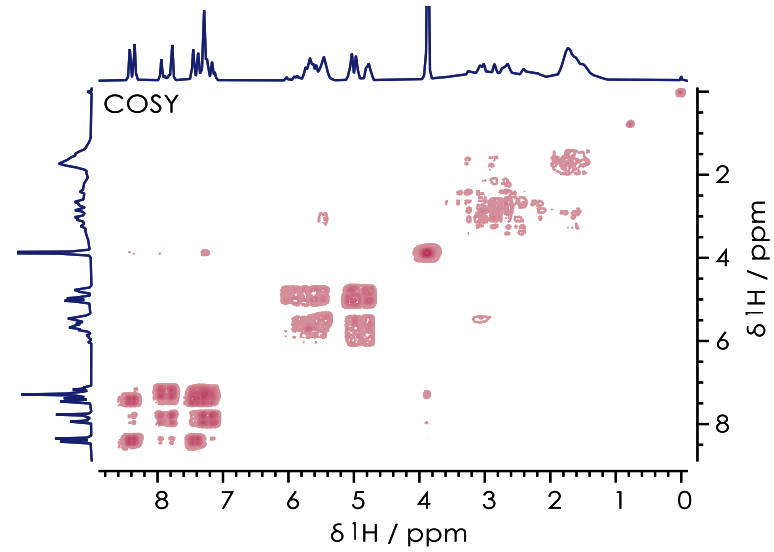

X-Pulse Spectra in NMRium



<https://nmr.oxinst.com/interactive-spectra-library>

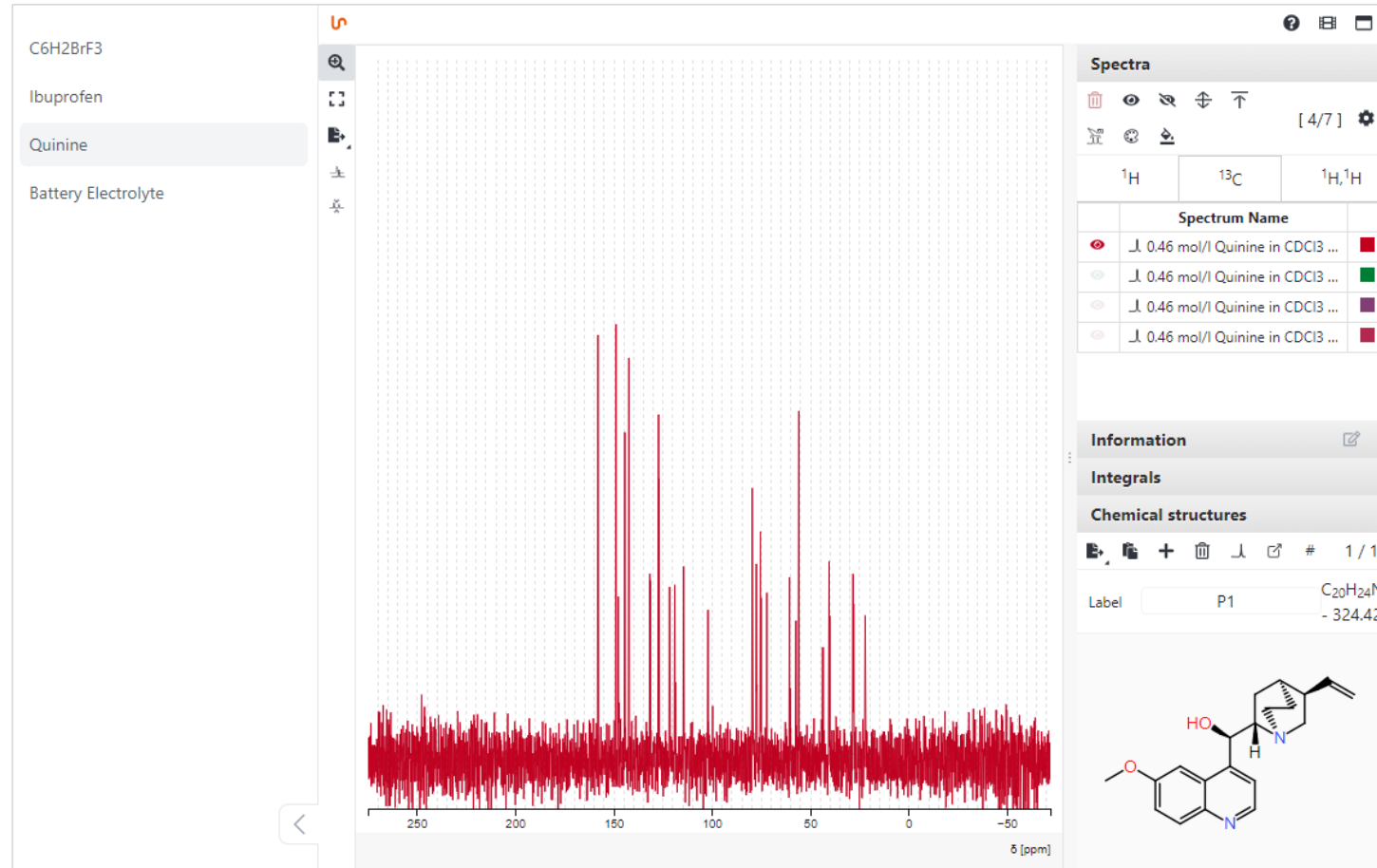


Reaction Monitoring & FlowNMR

www.oxinst.com/webinars/advances-in-reaction-monitoring-with-flownmr


Reaction Monitoring

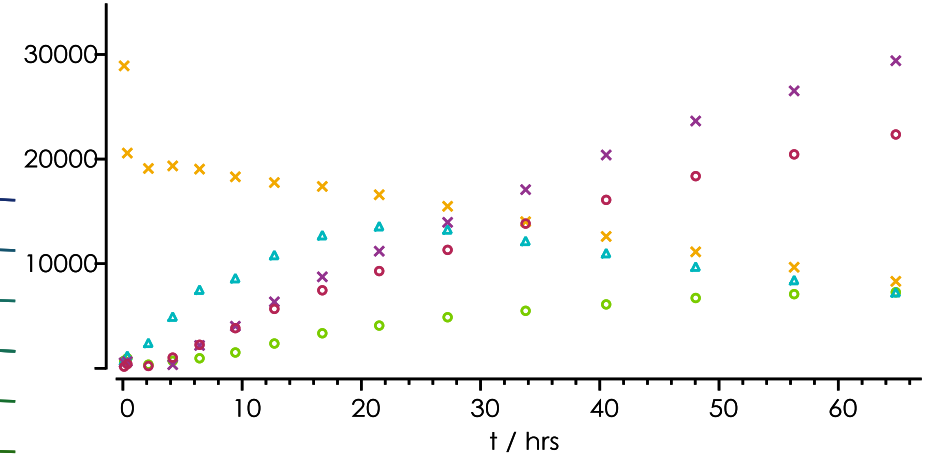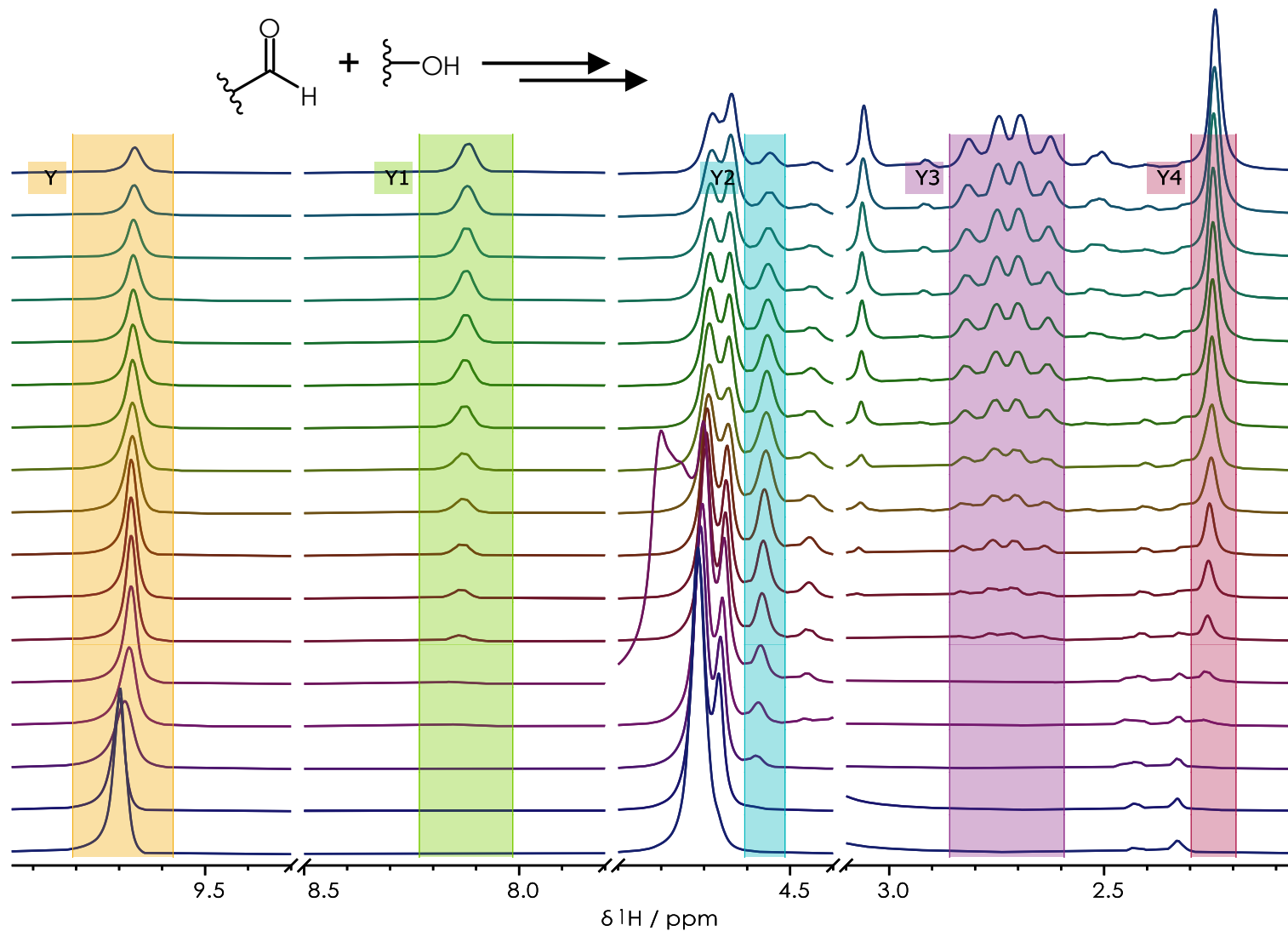

- The aldehyde signal at $\delta_H +9.7$ ppm, decreases as the starting material is consumed
- An intermediate species is observed at $\delta_H +4.6$ ppm
- Signals corresponding to product species can be identified, including singlets at $\delta_H +8.1$ and $+2.3$ ppm, and a quartet at $\delta_H +2.7$ ppm

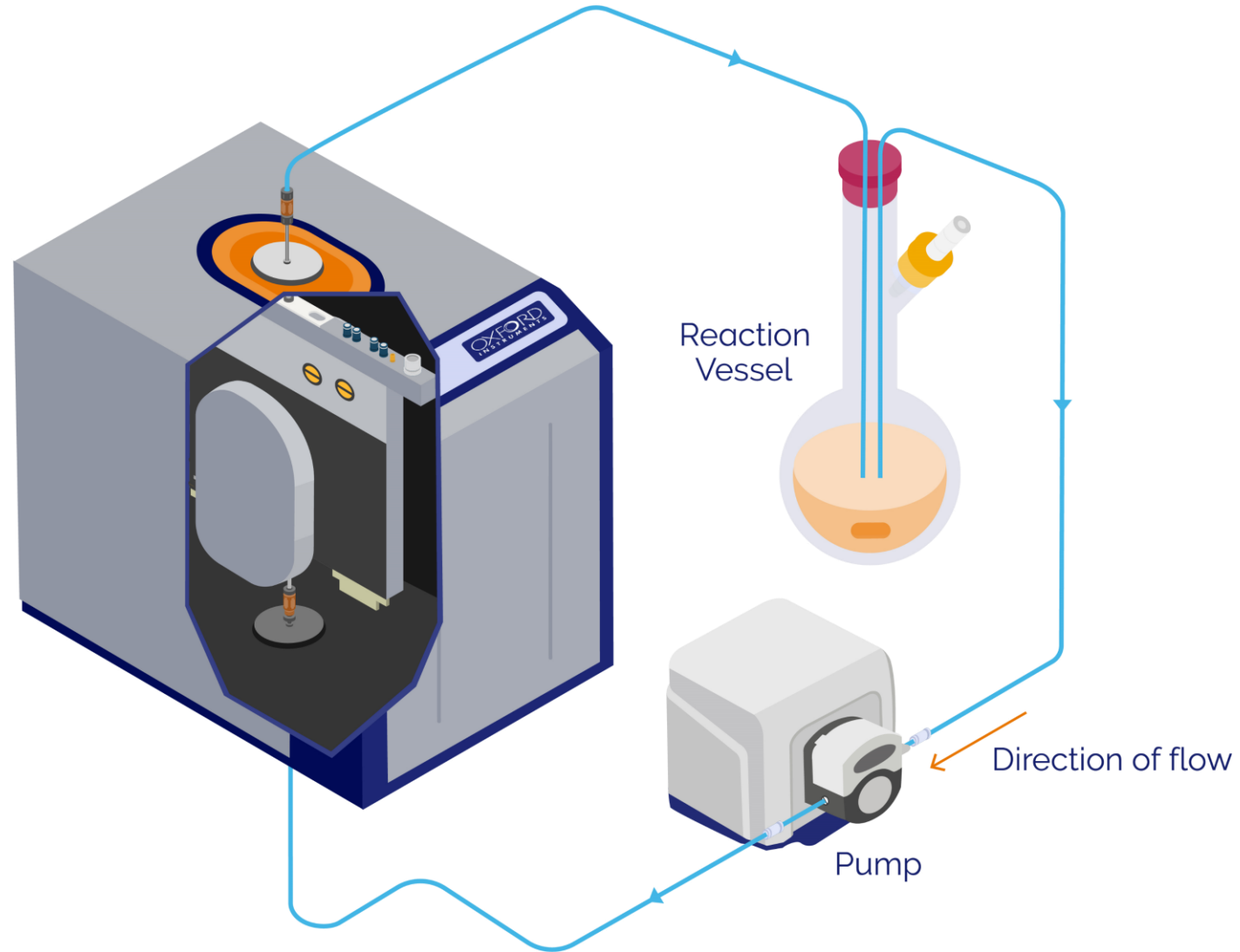


- Chemical reactions & processes can be monitored by flowNMR
 - calculation of rates
 - identification of intermediates / reaction completion
 - optimisation of reactions
- Wide range of applicable nuclei
 - ^1H , ^7Li , ^{11}B , ^{19}F , ^{23}Na , ^{31}P ...
- Quantitative NMR (qNMR) is possible in flow
 - quantification of $\text{Li}^+(\text{aq})$ and $\text{Na}^+(\text{aq})$ – extraction of lithium
- Three-axis field gradients allows for advanced experiments
 - solvent suppression (WET) in flow
 - measuring diffusion constants / polymer sizing in flow

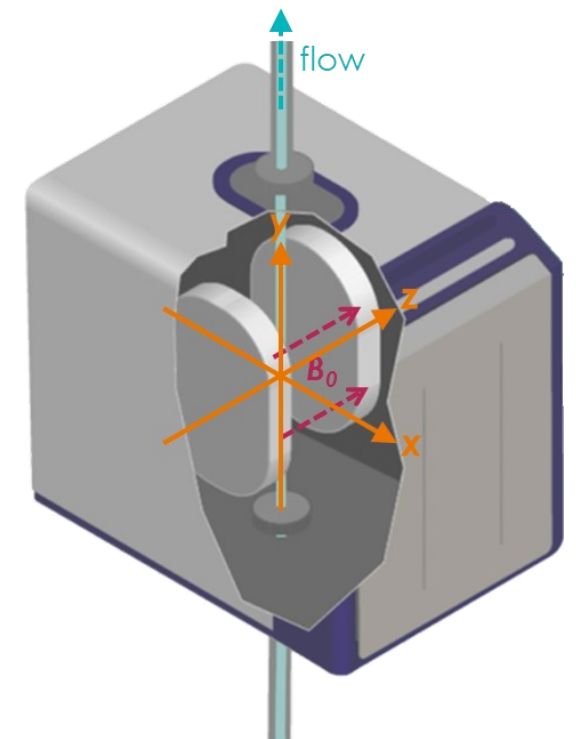

Time-Domain NMR

MQC+

What is Time-Domain NMR ?

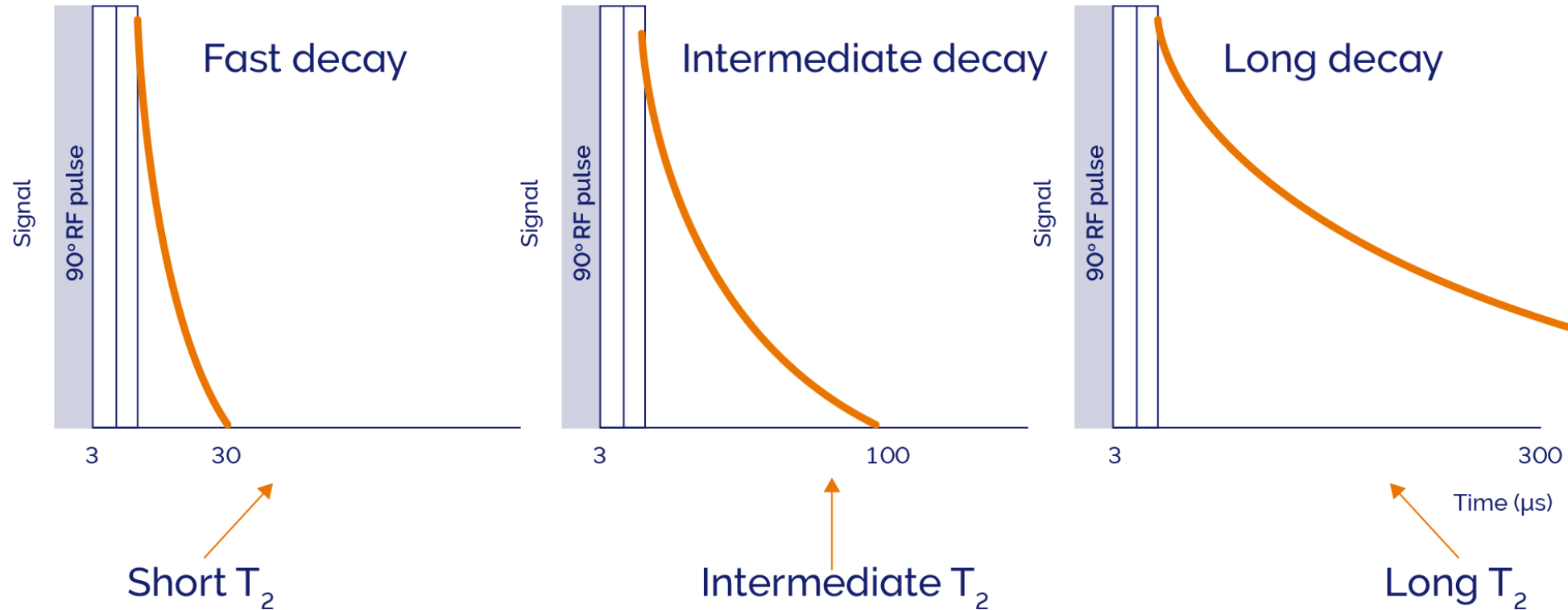

High Relaxation Rate

Low Relaxation Rate

Solids

- proteins, carbohydrates and fibre in foods
- polymer/biopolymer matrices of samples
- crystalline polymer components

Solid-like components

- bound water
- residual moisture
- amorphous polymer components

Liquids

- oils/fats
- spin-finish lubricants




- Oxford Instruments MQC family of TD-NMR Instruments
 - 5 MHz or 23 MHz permanent magnet
 - user removeable / exchangeable probes
 - Hydrogen-1 (proton) or Fluorine-19
 - 10 – 60 mm sample tube diameter
 - solid and/or liquid samples
 - *optional* variable temperature configuration (–5 to +70°C)
- QC/QA and full research grade systems
- Applications often replace ‘wet chemistry’ methods
- ISO & ASTM standard methods

Summary



- A good benchtop NMR spectrometer can (from an NMR point-of-view) do everything a high-field NMR spectrometer can do
 - that **doesn't** mean they can be used for all applications, though probably for more than you think ...
- Benchtop form-factor gives more system flexibility
 - fume cupboards, glove boxes, pilot plants ...

- NMR isn't just spectroscopy (and MRI) !
 - time-domain NMR / NMR relaxometry

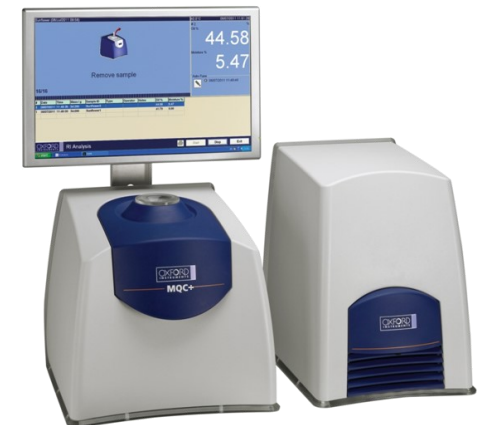

Any Questions ?

nmr.oxinst.com

www.linkedin.com/showcase/magnetic-resonance/

magres@oxinst.com

inclusive • innovative • trusted • purposeful